



www.labomed.com  
spectro@labomed.com

## 5MP Digital Camera for LB-356 Digital Zoom Stereo Microscope

### Easy to Install

Equipped with international standard C-type interface, it can be conveniently thread on both C-mount coupler and eyepiece adapter. It is compatible with the monocular, binocular, trinocular and any kind of microscopes, such as stereo, video, dissecting microscope or telescopes.

1. Connect video coupler to the device. Remove the dust cover and lens protection film, then connect the C-Mount coupler to the device.
2. Install the device on the microscope.



### Power Supply

Optional Plug Types: USA/ EU/UK and Japan.

Connect one end of DC adapter cable to the power port at the bottom of the device, and the adapter to 100-240V AC power outlet.

### Built-in Microscopy Software

After the device is turn on, touch the "YWCamera" Icon to open the camera application.

